

MICROLINKS TECHNOLOGY CO., LTD.

# Users Manual

---

**Installing MICROSCOPE Device Operates Manual**

**2019/01/19 Update Ver.3.0A**

## Contents

<b>1. Download &amp; Installation .....</b>	<b>3</b>
<b>2. Install Microscope application program.....</b>	<b>4</b>
<b>2.1 Install application program .....</b>	<b>4</b>
2.1.1 Start setup program.....	4
2.1.2 Welcome to the Microscope Setup Wizard.....	4
2.1.4 Now installing program .....	5
2.1.5 Complete install .....	6
2.1.6 After installation complete.....	6
<b>2.2 Uninstall MICROSCOPE AP .....</b>	<b>7</b>
2.2.1 Choose Microscope AP Uninstall see [Fig. 2-8].....	7
2.2.2 Confirm to Uninstall Microscope AP .....	7
2.2.3 Uninstalling Microscope AP .....	8
2.2.4 Finish Uninstall.....	8

## 1. Download & Installation

Please download the program from <http://www.vitiny.com> .

**Step 1: Uninstall MICROSCOPE AP. Please execute Autorun.exe, see Fig 1-1**



Fig 1-1

**Step 2: Install the program, see Fig 1-2 [ Refer to section 2.1]**

**Step 3: After complete installation, then press Exit.**



Fig.1-2 setup window

## 2. Install Microscope application program

Please select the image [Fig 1-1] to start setup application program.

### 2.1 Install application program

#### 2.1.1 Start setup program

Click setup application program, after it will pop up (please select a language) window on Fig.2-1, the initial language will follow up the pc OS system. If you want to cancel setup please click “Cancel” then click “OK”.

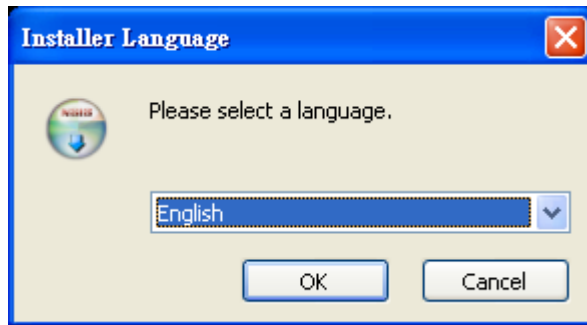


Fig. 2-1 setup language window

#### 2.1.2 Welcome to the Microscope Setup Wizard

After choosing the language, go to welcome to the MICROSCOPE setup wizard window see Fig.2-2, then please click “Next” then continue. If want to cancel please click “Cancel” button. In the folloing figs, “name” represents Microscope’s application program name, “X.XXX.XXXx” represents the application program version.

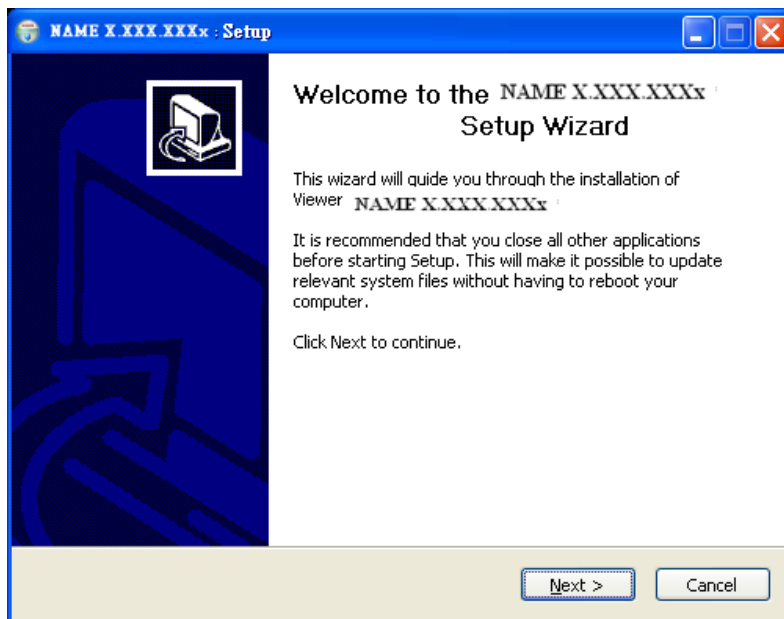


Fig. 2-2 welcome to the Microscope setup wizard

### 2.1.3 Choose installation route

Next step please choose the installation route window see [Fig.2-3]. If want to change file route click “Browse” to change setup path, then please click “install” then start installing.

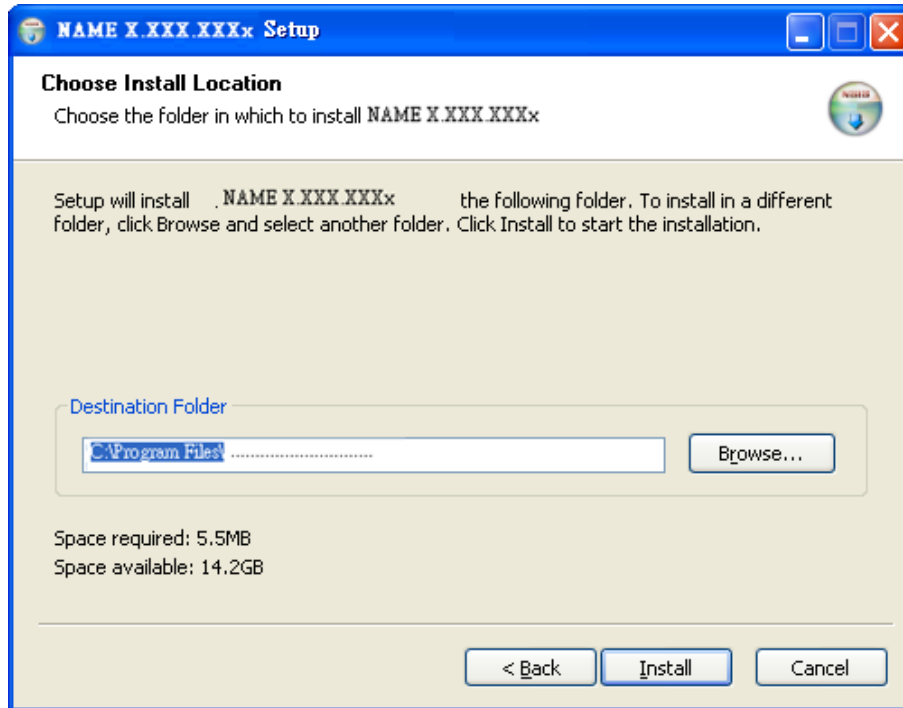


Fig. 2-3 choose installation route

### 2.1.4 Now installing program

Continue install wizard will start installing program as directory Fig..2-4.

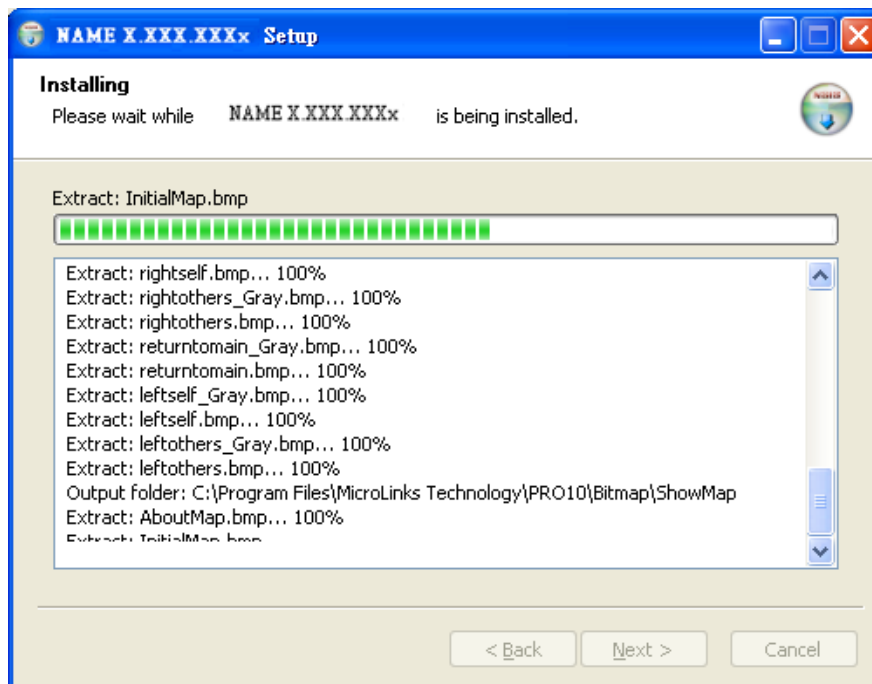


Fig.2-4 now installing

### 2.1.5 Complete install

When the program completes installation (Fig.2-5). Please click “Finish”.

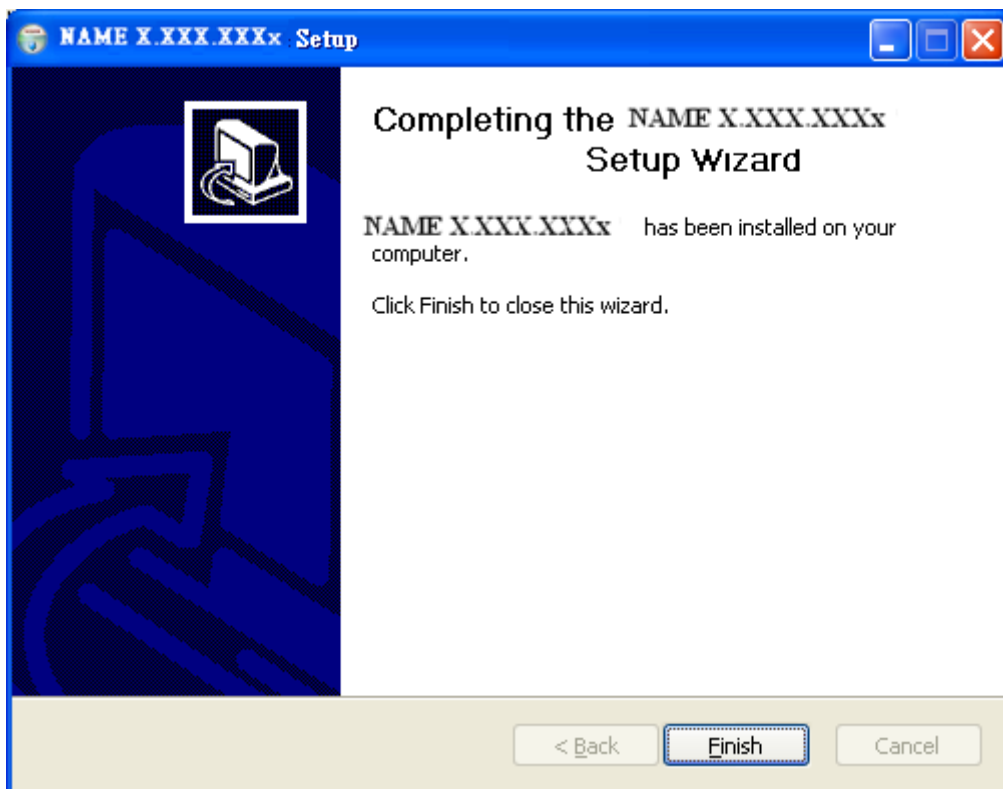


Fig.2-5 Microscope wizard install completed

### 2.1.6 After installation complete

Microscope application program shortcut will show on the Desktop as Fig.2-6.



Fig.2-6 Microscope shortcut

Or go “Start”-> All Programs-> MicroLinks Technology ->Microscope (see Fig 2-7).

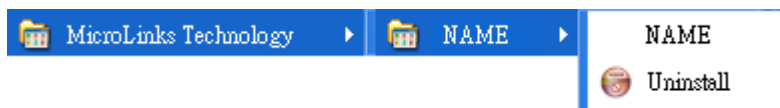


Fig. 2-7 The location of MICROSCOPE AP in all programs

## 2.2 Uninstall MICROSCOPE AP

2.2.1 Choose Microscope AP Uninstall see [Fig. 2-8]

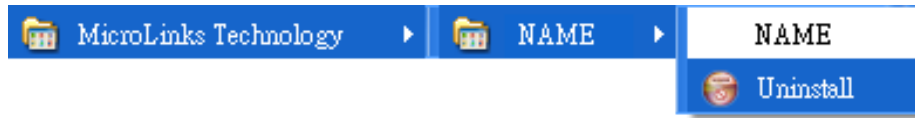


Fig. 2-8 Uninstall MICROSCOPE AP

2.2.2 Confirm to Uninstall Microscope AP

Please make sure if you want to completely remove Microscope application program.

See Fig 2-9.

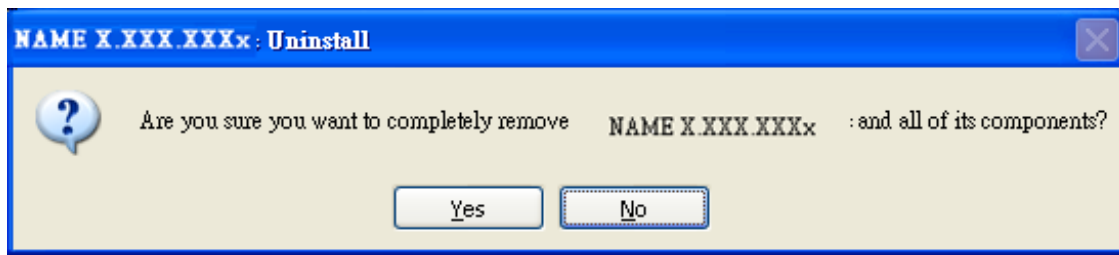


Fig. 2-9 Confirm uninstallation

### 2.2.3 Uninstalling Microscope AP

In processing of uninstall the Microscope AP, see Fig. 2-10

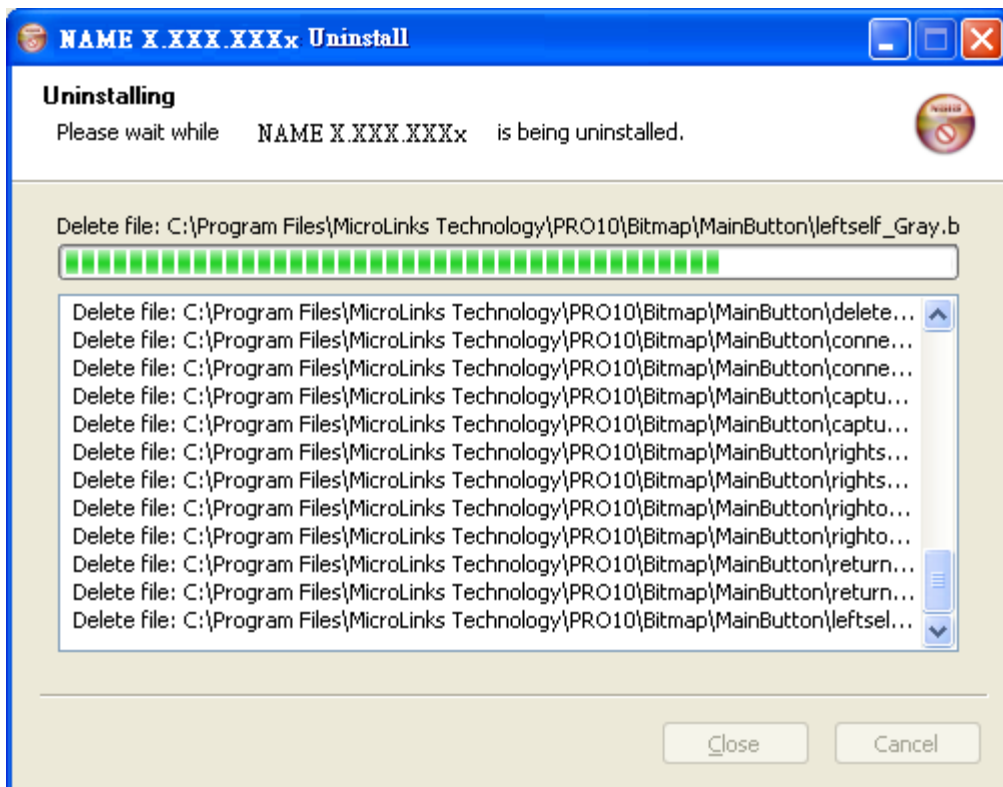


Fig.2-10 Uninstalling Microscope AP

### 2.2.4 Finish Uninstall

Fig. 2-11 shows that Microscope was successfully removed from your PC, please click OK to finish.



Fig. 2-11 Uninstall Completed